## INSTRUCTIONS

## **Biohazard Autoclave Gloves**

Catalog No. L1101



Gloves are suitable for temperatures up to 232°C (450°F). Always exercise caution when working with hot objects and equipment. Contact time with hot objects should always be minimized.

## DO NOT USE A WET GLOVE

## DO NOT HOLD GLOVE IN LIVE STEAM

Gloves will absorb liquids that conduct heat energy, which may cause burns to the hand.

Be careful when picking up sharp or pointed objects that could penetrate the glove.

Gloves are made of cotton terry cloth and can be machine washed and dried.

